



Product Information

SOLO+ Medium (SOL001)

Description

Selective enrichment media for use with Solus One *Listeria* assay.

Method for reconstitution

1. Dispense 59.7g of powder into 1L of distilled water.
2. Allow to soak for 10 minutes. Swirl to mix and dispense into suitable autoclavable containers.
3. Sterilize by autoclaving for 15 minutes at 121°C.
4. Allow to cool to ambient temperature.

Storage of Prepared Medium: Capped container 2 weeks at 2-8°C.

Sample enrichment

Warm SOLO+ to 30°C prior to use.

Solus Scientific Solutions Ltd

9 Mansfield Networkcentre, Millennium Business Park, Concorde Way, Mansfield, Nottinghamshire NG19 7JZ

Tel +44 (0)1623 429701 **Fax** +44 (0)1623 620977 **Email** enquiries@solusscientific.com

Web www.solusscientific.com

Registered Address: East Kilbride Laboratory, Torus Building, Rankine Avenue, East Kilbride, Glasgow, Lanarkshire G75 0QF Company Reg. No. SC361616, Registered in Scotland VAT No 982 2304 24