

# Palcam (PAC) Supplement

SUP011



For the isolation of *Listeria* spp.

## Presentation

This pack contains 10 single use vials containing sufficient supplement to give the correct working concentration when added to 500ml of Palcam broth (MED039).

## Shelf Life

This product has a shelf life of 2 years providing it is stored at 2-8°C. Shipping at ambient for up to 7 days is permissible. After rehydration some compounds or substances are unstable even when frozen and therefore unused supplement should be discarded.

## Contents

	Per vial	Conc. in medium
Polymyxin	5.0mg	10.0mg/L
Acriflavine	2.5mg	5.0mg/L
Ceftazidime	10.0mg	20.0mg/L

## Re-hydration

Wearing latex gloves aseptically reconstitute the contents of the vial with 5ml of sterile deionized water using a sterile pipette.

## Addition

Aseptically add the supplement to the media cooled to 47°C and immediately mix gently to evenly distribute.

## Safety Warnings and Precautions

If the vial is broken and the dry material or liquid escapes it, refer to appropriate SDS for possible hazards.